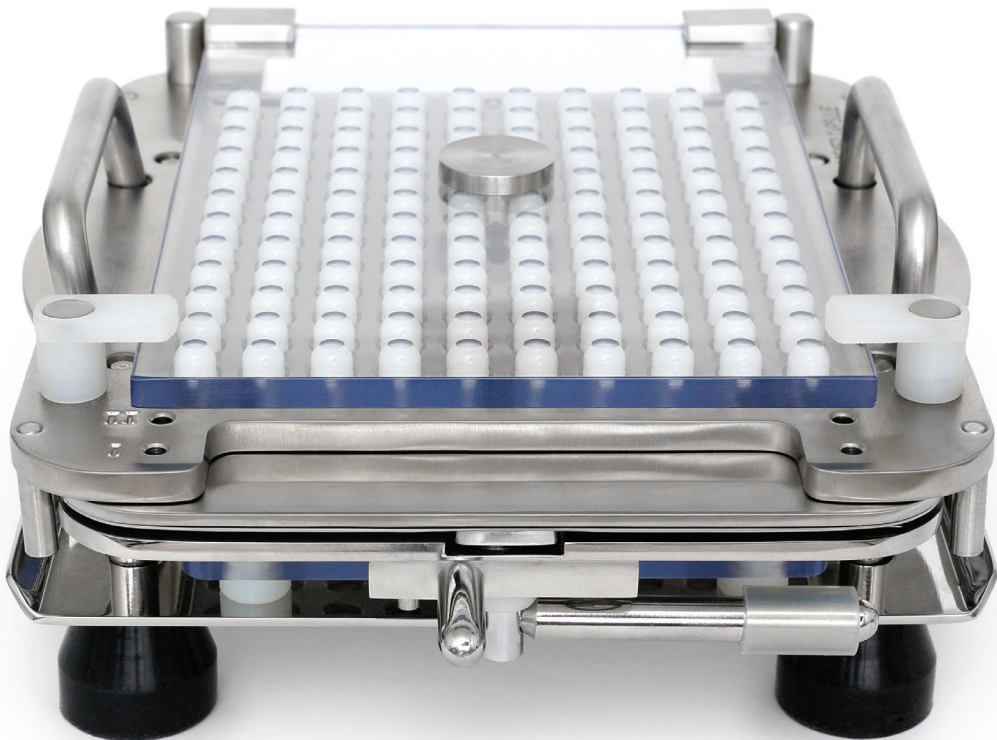
CapsiCards[®] System

Quick Start Manual
100 hole - 120 hole

V.1.0

Contents

Overview

Features: CapsiCards® System	5
Box Contents	10
Capsule Hole Number and Capsule Size Check	14

Initial Assembly / Capsule Size Change	16
---	-----------

Operating Instructions	18
-------------------------------------	-----------

CapsiCards® System Assortment	27
--	-----------

Materials and Cleaning

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CapsiCards®

Easy to use and flexible

With its innovative design, the CapsiCards® Machine is easy to handle and ensures quick, safe and hygienic capsule production.

Key features at a glance:

- Filler Base Kit available in 100 or 120 hole. Works with capsule sizes 00, 0, 1, 2, 3 and 4.
- Order Filler Change Parts for each capsule size when you need them.
- Inox sheets for reduced static, long-life and dishwasher cleaning.
- Adapter and Pusher speed capsule ejection from CapsiCards®.
- Fast lock cam design.
- Three plates for 100% capsule separation.
- Includes Powder Tray and Clamps to minimise powder loss.

- Includes Capsule Locker for fast and easy capsule locking.
- Size change requires no tools or adjustments.
- Easy cleaning as powder contact parts are Inox and dishwasher safe polycarbonate.
- Pharmacopoeia compliant for fill weight (USP/EP method).
- Filler Divider available for less than full batch of capsules.
- Powder Vibrator available for faster filling of hard-to-fill powders.
- System weights and dimensions available in the detailed Operating Manual (capsicardssystem.com and USB).

The CapsiCards® Machine has been developed especially for CapsiCards®.

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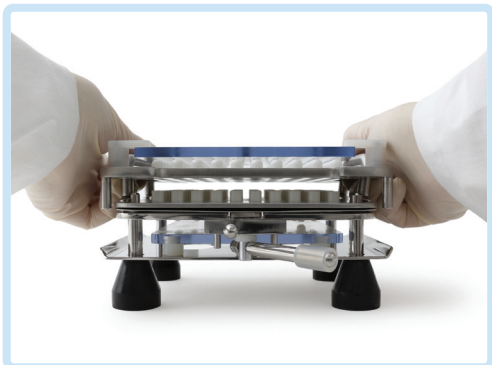
6 To separate capsules, press palms down on handles while lifting Caps Tray up with fingers. Remove Caps Tray with caps from Filler.

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7 Release Cam Lever so capsules fall down into filler.

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8 Place Powder Tray onto Filler. Optionally, attach Powder Tray Clamps.

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9 Pour powder onto Filler and spread. Use tapping, vibration or tamping to settle powder.

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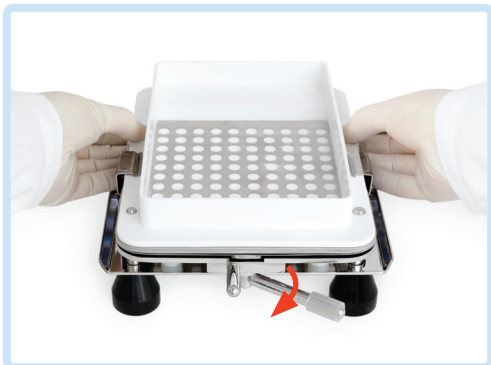
10a Tapping

Pull Cam Lever to gently hold bodies in place. Hold Filler Base Frame and gently tap against table a few times to settle powder. Powder Tray Clamps are recommended when tapping.

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10c Tamping

To fill additional powder, use Tamper.

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11 Remove Powder Tray. Place Caps Tray onto Filler.

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12 To lock capsules, press down on Locking Plate with thumbs while pulling up lifting plate. Repeat several times, moving hands from front to back along filler.

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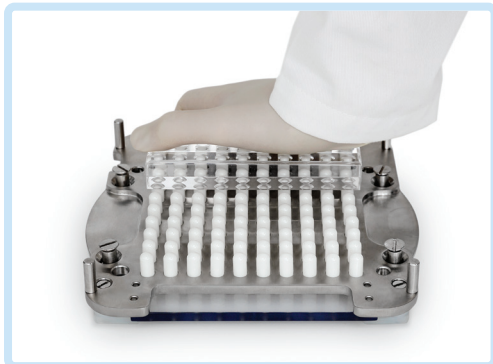
13 Remove Caps Tray and turn over onto table. Use Capsule Locker to lock capsules. Apply pressure until you feel capsules snap and lock.

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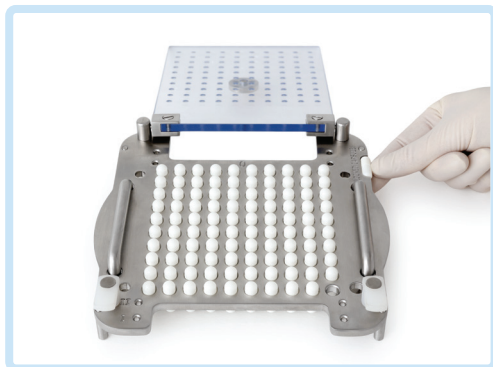
14 Capsules are fully locked. Check using locked capsule indicator.

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CapsiCards® System Assortment

CapsiCards® System

	Capsule no.	Capsule size	Material	Colour
CapsiCards® Machine – Filler Base Kit				
Filler Base Kit Filler Base, Powder Tray, Powder Tray Clamps, Powder Spreader, 5-pin Tamper, Capsule Locker, Pusher, Adapter	100 hole	Suitable for: 00, 0, 1, 2, 3, 4	Inox (Stainless Steel 316)	n/a
Filler Base Kit Filler Base, Powder Tray, Powder Tray Clamps, Powder Spreader, 5-pin Tamper, Capsule Locker, Pusher, Adapter	120 hole	Suitable for: 00, 0, 1, 2, 3, 4	Inox (Stainless Steel 316)	n/a
CapsiCards® Machine – Filler Change Parts				
Filler Change Parts Body Sheets (3), Caps Tray with Locking Plate, Spacer Plate	100 hole	Available in: 00, 0, 1, 2, 3, 4	Inox (Stainless Steel 316)	n/a
Filler Change Parts Body Sheets (3), Caps Tray with Locking Plate, Spacer Plate	120 hole	Available in: 00, 0, 1, 2, 3, 4	Inox (Stainless Steel 316)	n/a
CapsiCards®				
CapsiCards®	50 capsules/card	Available in: 00, 0, 1, 2, 3, 4	Hard gelatin	Clear, white, ivory
CapsiCards®	60 capsules/card	Available in: 00, 0, 1, 2, 3, 4	Hard gelatin	Clear, white, ivory

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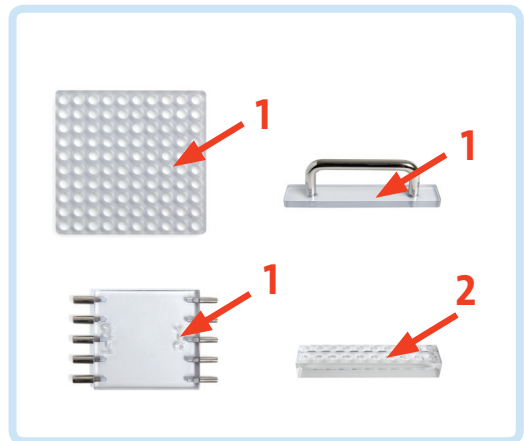
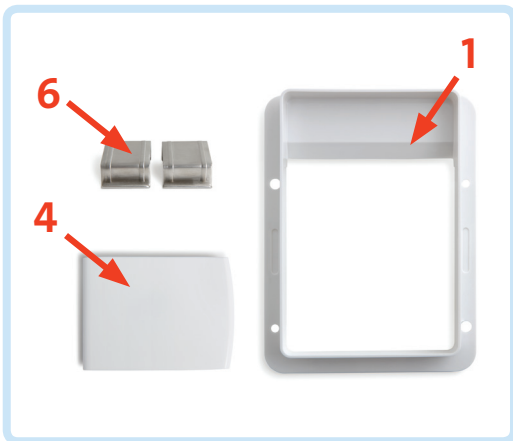
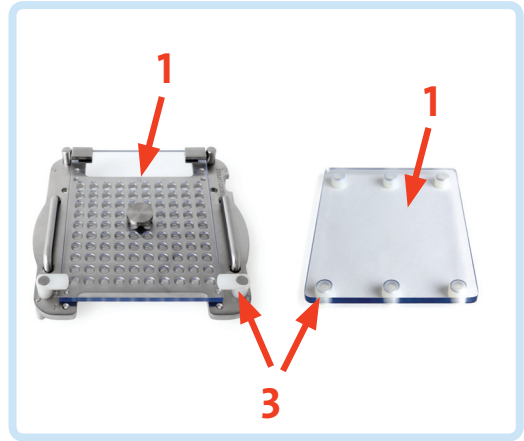
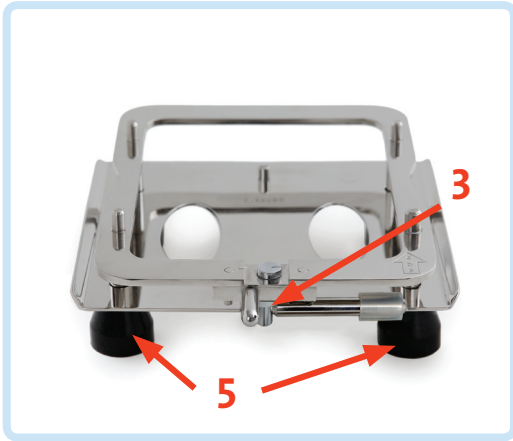
Cleaning

Automatic dishwasher

All parts are fully dishwasher safe. Plastic parts should be placed in top rack of dishwasher (< 70 °C / 160 °F).

Cleaning agents

We recommend cleaning using household hand dishwashing liquid. All items may be cleaned with NaOH (caustic soda) 1-2% solution, 1% sodium thiosulphate solution, or liquid detergents with up to 4% sodium hypochlorite.

Isopropyl alcohol

Wipe parts with a cloth dipped in isopropyl alcohol, then immediately wipe dry with a clean cloth. Do not use on acrylic capsule locker. Acrylic parts can be cleaned with a solution of 15% ethyl alcohol (ethanol) and 85% water.

Hot air drying

Up to 60 °C / 140 °F. Remove plastic parts from oven after 15-20 minutes. If using portable hair dryer, keep a distance of 30 cm (12 inches) from all items.

Autoclave

Do not autoclave capsule locker or powder spreader. Other parts may be autoclaved at 121 °C / 250 °F, 15 psi pressure for 15 minutes.

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Optional Accessories

Partial Batch Filler Divider

Fill fewer than the maximum capsules per batch.
Use Filler Divider to block empty holes during powder filling.
Works with 100 and 120 hole fillers.

Material:

Inox (SS316), PET and silicone rubber.

How to use:

1. Cut CapsiCards® to desired number of capsules.
2. Follow steps 1 - 7. See Chapter Operating Instructions, page 18.
3. Place Powder Tray onto Filler.
4. Place Filler Divider onto Filler to cover empty holes.
5. Follow steps 9 - 14. See Chapter Operating Instructions, page 21.

Vibrator

Recommended for faster filling of powders that bridge or are granular, fluid or flour-like. Reduces weight variation of fine powders. Works with 100 and 120 hole fillers.

100-120 Volts / 50-60Hz or 220-240 Volts / 50Hz.

How to use:

Place Filler on Vibrator. Use vibration to pack powder in step 10. See Chapter Operating Instructions, page 23.

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User Tips

More details available in detailed Operating Manual (capsicardssystem.com and USB).

Capsule bodies not level with body sheet set after separation

It is normal for the top of the capsule bodies to sit unevenly in the filler after separation from the caps. Capsules have slight variation in diameter and height, which allows some to fall through to the bottom plate while others are held by the body sheets. No capsules should sit higher than the top sheet. If they do, use your hand to pat them into place before pouring powder. After spreading powder and doing your first tapping, vibration, or tamping of the powder, the capsule bodies will settle properly into place and sit 0.5 to 1 mm lower than the top surface of the body sheet set.

Capsules are difficult to lock

Depending on the powder, up to 1 kg of force is required to fully lock one capsule. For 120 hole machines, the force required to lock all 120 capsules simultaneously could be as much as 100 kg. If, due to operator strength or powder properties, enough force cannot be achieved, use of the capsule locker will ensure fully locked capsules.

Capsules are too soft or too brittle

Capsule storage and filling should be from 15 to 25 °C (59 to 72 °F) with relative humidity from 35 to 65%.

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Declaration of Conformity

DECLARATION OF CONFORMITY

Product name Hand operated capsule filling machine
Model CapsiCards® Machine
Manufacturer Custom Capsules Pvt. Ltd.
B 8/1 Camlin Naka, Tarapur MIDC, Boisar,
Dist: Thane, Maharashtra, India, 401506

Contact person **Customer support** CE certification questions
For contact details of Mr. A.H. Varma
your local supplier, Phone: +91 2525-272538
please see page 46. Fax: +91 2525-272537

It is hereby confirmed that the above product complies with the requirements set out in the following directive:

2006/42/EC Machinery directive

The harmonized standards applied are as follows:

EN292/1 & EN292/2:1991 Safety of machinery - Basic concepts and general principles for design.

EN349:1993 Safety of machinery - Minimum gaps to avoid crushing of parts of the human body.

I hereby declare that the above product has been designed to comply with the relevant sections of the above referenced standards. The unit complies with all applicable essential requirements of the directive.

The product carries the CE mark.

Signed: 

Place: MIDC Tarapur, India

Name: A.H. Varma
Position: Assistant manager Q.A.

Date: 14 Jan. 2014

CapsiCards® is a trademark of Fagron Group BV.

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Conformiteitsverklaring

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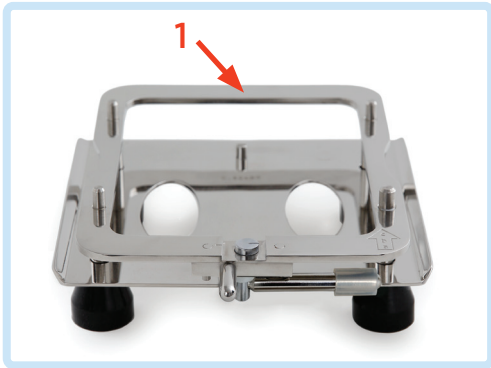
Warranty

12-month parts replacement in case of manufacturing defects. On-site service is not included.

The warranty does not cover damage caused by accident, misuse, alterations, tampering, negligence, or operation that does not conform to the instructions in the operating manual.

Please include Filler Base Serial Number with all correspondence (1).

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Manufacturer: Custom Capsules Pvt. Ltd. for Torpac Inc. USA. Made in India.