


Item Code: P1X PBK 00-4 for sizes 00-4 & 0e1-2e1







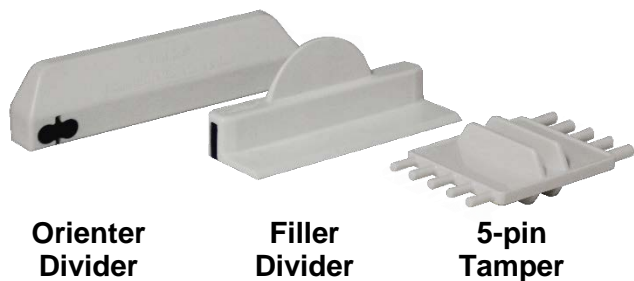

**Orienter
Divider**

**Filler
Divider**

**5-pin
Tamper**

Parts shipped are of new and improved design (see above).
Parts shown in instructions below are of original design.

Easiest way to fill less than a full batch!

-  Orienter Divider and Filler Divider design makes the PBK very easy to use. Takes seconds to install and remove.
-  Eliminates need for tape, cardboard and other non-GMP methods for making less than a full batch.
-  Works with ProFiller 1000 series standard white polycarbonate Powder Tray.
-  Prepare batch sizes in increments of 10 capsules.
-  Made of polycarbonate: dishwasher safe, isopropyl alcohol safe.
-  Autoclave safe at 121°C, 15 psi pressure for 30 Minutes.


Instructions for Use

Batch sizes

ProFiller 1060: 10-30 capsules



ProFiller 1100: 10-70 capsules

ProFiller 1120: 10-90 capsules

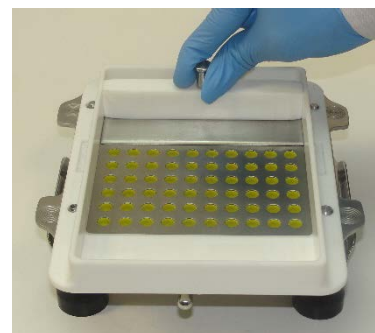

1. Place Orienter Divider on Orienter at half the desired batch size¹. Orient capsules in positions I and II as usual.

2. Separate capsules in Filler as usual. Place Powder Tray onto Filler. Place Filler Divider onto Filler.

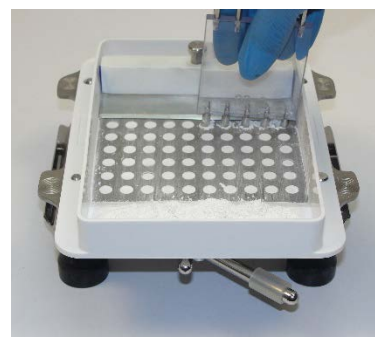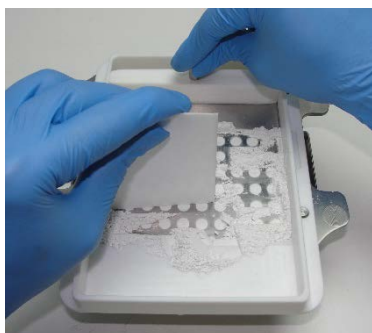

3. Pour powder and spread using Powder Spreader. If needed, press down on Divider to reduce powder loss.

4. If using a 5-Pin Tamper or tapping, gently pull Cam Lever toward post to secure bodies in Filler.

5. Use regular Tamper or 5-pin Tamper to tamp powder, or tap or vibrate as usual. Release Cam Lever. Remove Powder Tray and lock capsules as usual.

Other Accessories



Tap Densimeter



Powder Vibrator

For more information visit
www.profiller.com

¹ Orienter Divider works for all sizes regardless of Orienter Tray ramp direction.